# LETTER of ENDOTOXIN

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### Low Endotoxin Recovery (LER) について(1)

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1. Chen, J. Low Endotoxin Recovery in Common Biologics Products. Presented at PDA 8<sup>th</sup> Annual Global Conference on Pharmaceutical Microbiology, Bethesda, MD, (2013).
2. 高岡文ら、微量金属イオンがエンドトキシンのリムルス活性に及ぼす影響、日本薬学会 110 年会、(1990).
3. Patricia, F. Endotoxin Challenge – A Regulatory Perspective. Presented at PDA 9<sup>th</sup> Annual Global Conference on Pharmaceutical Microbiology, Las Vegas, NM, (2014)
4. Cooper, J.F. How Can Water Interference with BET?. Presented at the Pharmaceutical Microbiology Forum Bacterial Endotoxins Summit Meeting, Philadelphia, PA, (2014).
5. Platco, C. Lab Experience Low Endotoxin Recovery. Presented at the Pharmaceutical Microbiology Forum

Bacterial Endotoxins Summit Meeting, Philadelphia, PA, (2014).

6. Bolden, J. et al. Evidence Against a Bacterial Endotoxin Masking Effect in Biologic Drug Product by Limulus Amebocyte Lysate Detection. *PDA J Pharm Sci and Tech.* **68**, 472-447, (2014)
7. Reich, J. Heterogeneity of Potential Endotoxin Contaminations in Parenteral Drugs. Presented at the Parenteral Drug Association Conference, Berlin, Germany, (2015).
8. Tsuchiya, M. Possible Mechanism of Low Endotoxin Recovery. *American Pharmaceutical Review.* **17**, 1-5.(2014)
9. Reich, J. New Control Strategies for Monitoring of Endotoxin Masking in Biologics. Presented at the Parenteral Drug Association Conference, Berlin, Germany, (2014).